

# NOTICE

**When opening the autoclave the following PPE is required:**

- Eye protection (Safety glasses or goggles)
- Lab Coat
- Thermal Gloves
- Use the door as a shield when opening.

**The following should not go in the autoclave:**

- Aerosols
- Chemicals (corrosives, flammables, volatiles, cleaning products)
- Radioactive materials

# NOTICE

**When opening the autoclave the following PPE is required:**

- Eye protection (Safety glasses or goggles)
- Lab Coat
- Thermal Gloves
- Use the door as a shield when opening.

**The following should not go in the autoclave:**

- Aerosols
- Chemicals (corrosives, flammables, volatiles, cleaning products)
- Radioactive materials